



Micropipette User Guide

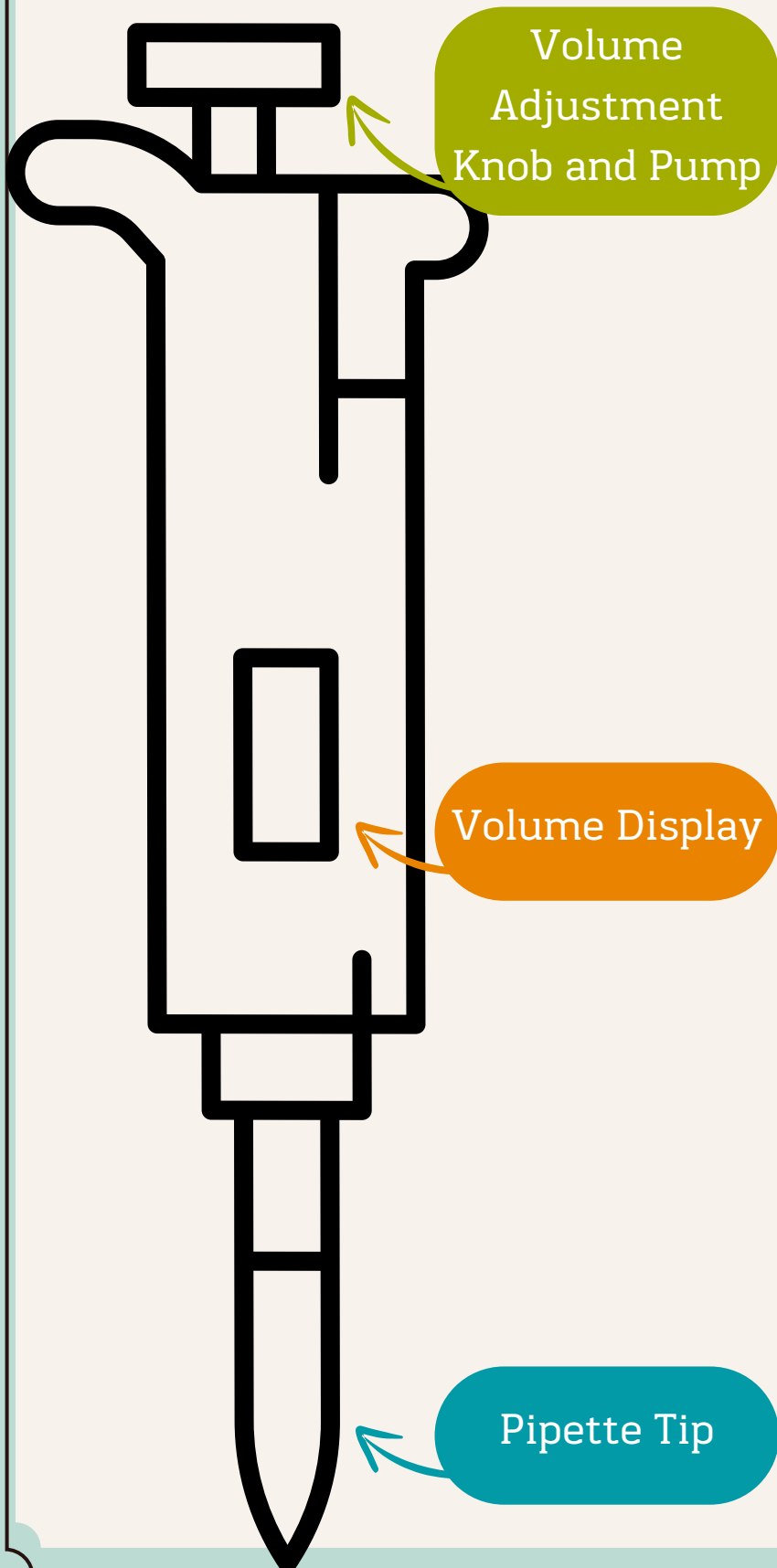


1. Insert the micropipette into the tip and press firmly. Twist back and forth a couple of times to make sure it's fitted snugly.
2. Turn the volume adjustment knob until the volume display shows the amount you want to dispense. Turn clockwise to decrease the volume and counter-clockwise to increase the volume.
3. Insert the tip of the pipette into the liquid you want to draw up.
4. Hold the pipette vertically, not at an angle, while drawing up liquid using the chart on the other side of this sheet.


Micropipette User Guide

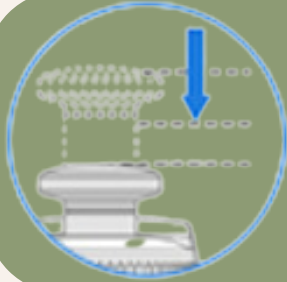

Before inserting the micropipette into the liquid, press the pump down to the first stopping point.

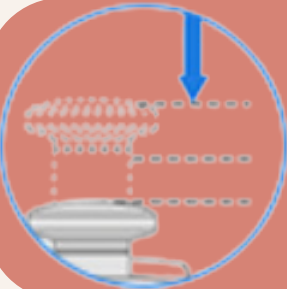

Dip the pipette tip into the liquid and slowly release to the original point.

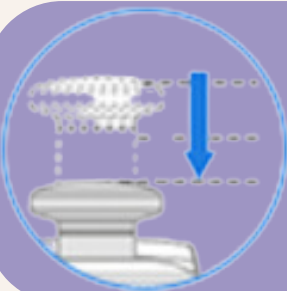

Dispense the liquid by pressing the pump in to the second stop. Hold until all liquid is dispensed.